



Table of Contents

Goals of this Experimental Guide	3
What is Patch-Clamp Electrophysiology and Why is it Important?	3
Who This Guide Can Help	5
Contact Precisionary Instruments for Any Additional Experimental Support	6
Gross Anatomic Location of Mouse Brain Structures	7
Mouse Brain Anatomy	7
Key Differences and Similarities	7
Human Brain Anatomy	9
Using Human Brain Tissue in Patch-Clamp Studies	10
Essential Materials and Reagents for Patch-Clamp Electrophysiology	11
Solutions & Recipes for Patch-Clamp Electrophysiology	12
External Solutions	12
Internal Solutions	13
Final Tips and Hints for Making External and Internal Solutions for Patch-Clamping	15
Importance of Osmolarity	15
Practical Solution Preparation	15
Additional Tips	
Helpful Hints	16
Agarose Solution	17
70% Ethanol	18
Making Acute Brain Slices from Mouse	18
Preparing for Slicing Mouse Brain Slices	18
Dissection of the Mouse Brain for Slicing	20
Sectioning Acute Brain Slices on the Compressstome Vibratome	21
Embedding Brain Tissue into a Specimen Tube: A Step-by-Step Guide	21
Cutting Acute Brain Slices on the Compressstome Vibratome	23
Collecting and Incubating Acute Brain Slices	24
Recording from Acute Brain Slices	25
Making Recording Pipettes	26
Lowering the Pipette Towards the Target Cell	26
Making the Gigaseal	28
Establishing the Type of Patch-Clamp Recording	29
Whole-Cell Recording	29
Perforated Patch Recording	30
Single Ion Channel Recordings	32
What to Include in Your Ephys Logbook	33
Making Bipolar Electrodes for Stimulation	34
Troubleshooting Patch-Clamp Electrophysiology	37
More Questions? Contact Us!	39

